12 RESPONSIBLE
CONSUMPTION
AND PRODUCTION

Vestfrost Solutions is working towards reaching the UN - Global Sustainable Development Goals by 2030.

The Sustainable Development Goals are the blueprint to achieve a better and more sustainable future for all. In order to implement Goal no 12 "Responsible Consumption and Production", this manual has been printed on recycled paper.

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WARNING

As the appliance contains flammable refrigerant, as stated on nameplate, it is essential to ensure that the refrigerant pipes are not damaged.

The quantity and type of the refrigerant used in your appliance is indicated on the rating plate.

Standard EN378 specifies that the room in which you install your appliance must have a volume of 1m³ per 8 g of hydrocarbon refrigerant used in the appliances. This is to avoid the formation of flammable gas/air mixtures in the room where the appliance is located in the event of a leak in the refrigerant circuit.

WARNING:

Keep ventilation openings in the appliance or in built-in structures must be kept clear.

WARNING:

Do not use other mechanical devices or means to accelerate the defrosting process or to remove rime other than those recommended by the manufacturer.

WARNING:

Do not damage the refrigerant system.

WARNING:

The appliance may not be exposed to rain.

WARNING:

This appliance is not intended for use by persons (including children) with reduced physical, sensory or mental capabilities, or lack experience and knowledge, unless they have been given supervision or instructions concerning use of the appliance by a person responsible for their safety. Children should be supervised to ensure that they do not play with the appliance.

WARNING:

Children must not play with, on, or around the appliance.

WARNING:

Children must not clean the appliance or carry out general maintenance unless they are at least 8 years old and are being supervised.

WARNING:

Always, keep the keys in a separate place and out of reach of children.

WARNING:

Before servicing or cleaning the appliance, switch off circuit breaker.

**WARNING:**

Danger risk of fire or explosion. Flammable refrigerant used, as stated on nameplate. To be repaired only by trained personnel.

WARNING:

Do not store explosive substances such as aerosol cans with a flammable propellant in this appliance.

WARNING:

When positioning the appliance, ensure the power cord is not trapped or damaged.

WARNING:

Do not locate multiple portable socket-outlets or portable power supplies at the rear of the appliance.

WARNING:

Appliance use flammable insulation blowing gas. For information about safe disposal, please contact your local disposal service. See section for Disposal.

Contents

WARNING	2	Starting device replacement for VLS504A AC.....	32
Periodic preventive maintenance checks ...	5	Run capacitor replacement VLS204 to 404A AC.....	35
Complete spare part list VLS204A AC	6	Run capacitor replacement for VLS504A AC.....	36
Complete spare part list VLS304A AC	8	Thermometer replacement	37
Complete spare part list VLS354A AC	10	Fan replacement VLS504A AC	40
Complete spare part list VLS404A AC	12	Compressor replacement	41
Complete spare part list VLS504A AC	14	On-site checklist	42
Vital components	16	Trouble shooting	43
Health and safety guidance – Warning!.....	17	Technical support.....	44
Required basic tools	18	Recycling procedures	45
Wire diagram VLS204A AC / 304A AC	19		
Wire diagram VLS354A AC / 404A AC	20		
Wire diagram VLS504A AC	21		
Motor compartment	22		
Thermostat replacement.....	23		
Thermostat adjustment.....	26		
Thermostat sensor replacement.....	27		

PQS Code	Model	PQS Performance specifications Specification reference:	PQS Independent type-testing protocol Product verification protocol:
E003/109	VLS 204A AC	E003/RF03.4	E003/RF03-VP.3
E003/110	VLS 304A AC	E003/RF03.4	E003/RF03-VP.3
E003/111	VLS 354A AC	E003/RF03.4	E003/RF03-VP.3
E003/112	VLS 404A AC	E003/RF03.4	E003/RF03-VP.3
E003/113	VLS 504A AC	E003/RF03.4	E003/RF03-VP.3

Periodic preventive maintenance checks

Daily Check:

Monitor Temperature.
Lid fits and lock tight to cabinet
Lid gasket not faulty.

Weekly maintenance:

Remove any water at the bottom of the refrigerator with a cloth.
Wipe of water droplets on the inside wall.

Monthly maintenance:

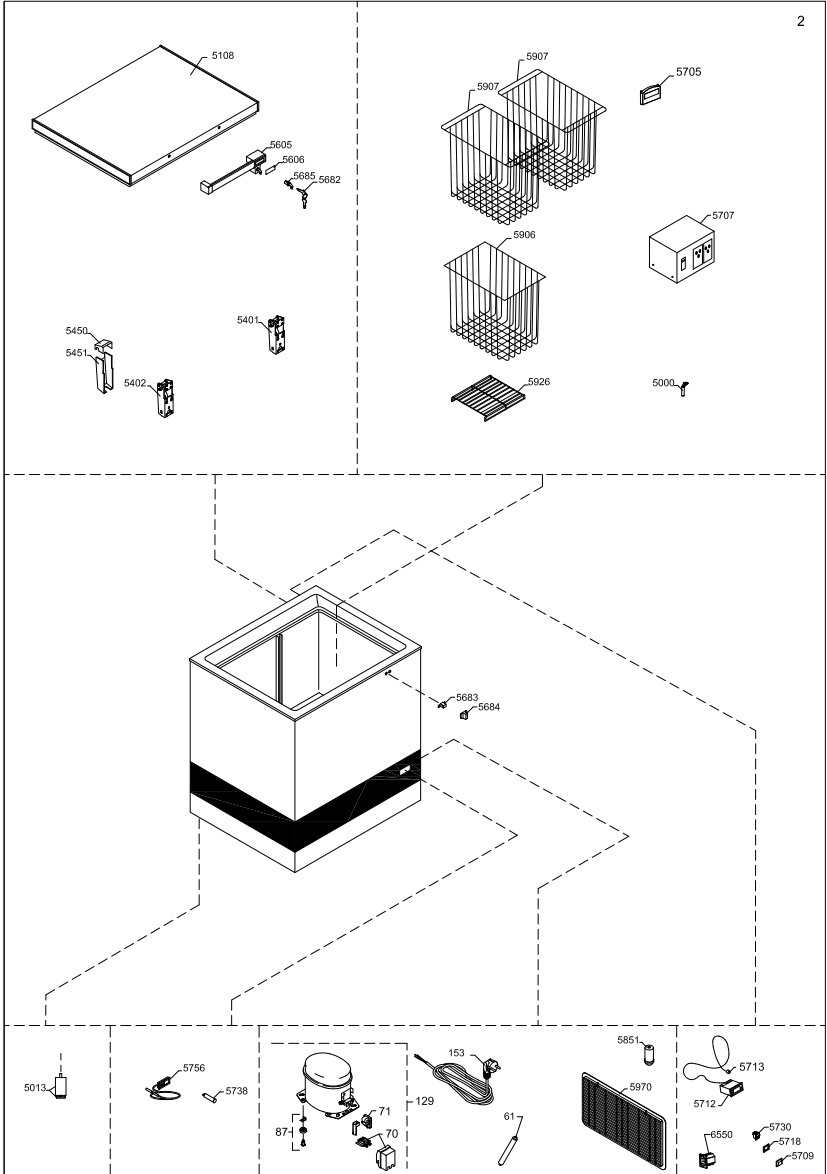
Clean grille for compressor compartment.
Clean the refrigerator with lukewarm water and mild detergent.

Yearly maintenance:

Check electrical connections and components.


Complete spare part list VLS204A AC

Position	Item number	Item name
0061	0-6538001	Filter drier,
0070	A921001	Cover + wiring clamp
0087	0-6038175	Base plate fittings, complete
0129	8-036038297	Compressor - TLES7KTK , Complete,
0153	7535254	Power cord EU H05VV-F 3G1 C13 jack 90°
5000	3010049	Drain plug
5013	3040400	Adjustable foot
5108	5000917680	Lid foamed without handle/hinges
5401	1510133	Hinge
5402	1510136	Hinge with spring
5450	3011135-01	Top part for hinge cover
5451	3010032-01	Bottom part for hinge cover
5605	304090501	Handle with lock (push and turn)
5606	8090342-94	Inlay for handle,
5682	1510238	Nikel-plated Key
5683	2040145	Catch for handle
5684	3010265-01	Cover for catch
5685	600098801	Lock with keys,(push and turn)
5705	7020406	Temperature monitoring device Fridge-Tag
5707	7020248	Voltage stabilizer
5709	7060104	Frame for cover rocker switch bezel
5712	7095514	Electronic controller ERC 111A
5713	7020960	Sensor NTC S1, Cool
5718	7060105	Rocker Switch Cover
5730	7020245	"on-off" switch
5738	7010138	Pilot lamp, green
5756	7020392-03	Thermometer, solar
5810	A921115	Starting device, 103N0021
5851	6520229	Run capacitor
5906	3510016	Basket 275 mm
5907	3510015	Basket 236 mm
5926	3510032	Bottom grating
5970	3010308-01	Motor screen
6550	7080143	Connector




Complete spare part list VLS304A AC

<u>Position</u>	<u>Item number</u>	<u>Item name</u>
0061	0-6538001	Filter drier,
0070	A921001	Cover + wirring clamp
0087	0-6038175	Base plate fittings, complete
0129	8-036038297	Compressor - TLES7KTK , Complete,
0153	7535254	Power cord EU H05VV-F 3G1 C13 jack 90°
5000	3010049	Drain plug
5013	3040400	Adjustable foot
5112	50009112163	Lid foamed without handle/hinges
5401	1510133	Hinge
5402	1510137	Hinge with spring
5450	3011135-01	Top part for hinge cover
5451	3010032-01	Bottom part for hinge cover
5605	304090501	Handle with lock (push and turn)
5606	8090342-94	Inlay for handle,
5682	1510238	Nikel-plated Key
5683	2040145	Catch for handle
5684	3010265-01	Cover for catch
5685	600098801	Lock with keys,(push and turn)
5705	7020406	Temperature monitoring device Fridge-Tag :
5707	7020248	Voltage stabilizer
5709	7060104	Frame for cover rocker switch bezel
5712	7095514	Electronic controller ERC 111A
5713	7020960	Sensor NTC S1, Cool
5718	7060105	Rocker Switch Cover
5730	7020245	"on-off" switch
5738	7010138	Pilot lamp, green
5756	7020392-03	Thermometer, solar
5810	A921115	Starting device, 103N0021
5851	6520229	Run capacitor
5874	A921004	Cover + wirring clamp
5906	3510016	Basket 275 mm
5907	3510015	Basket 236 mm
5919	3520056	Basket 225 mm
5925	3510031	Bottom grating
5970	3010308-01	Motor screen
6550	7080143	Connector

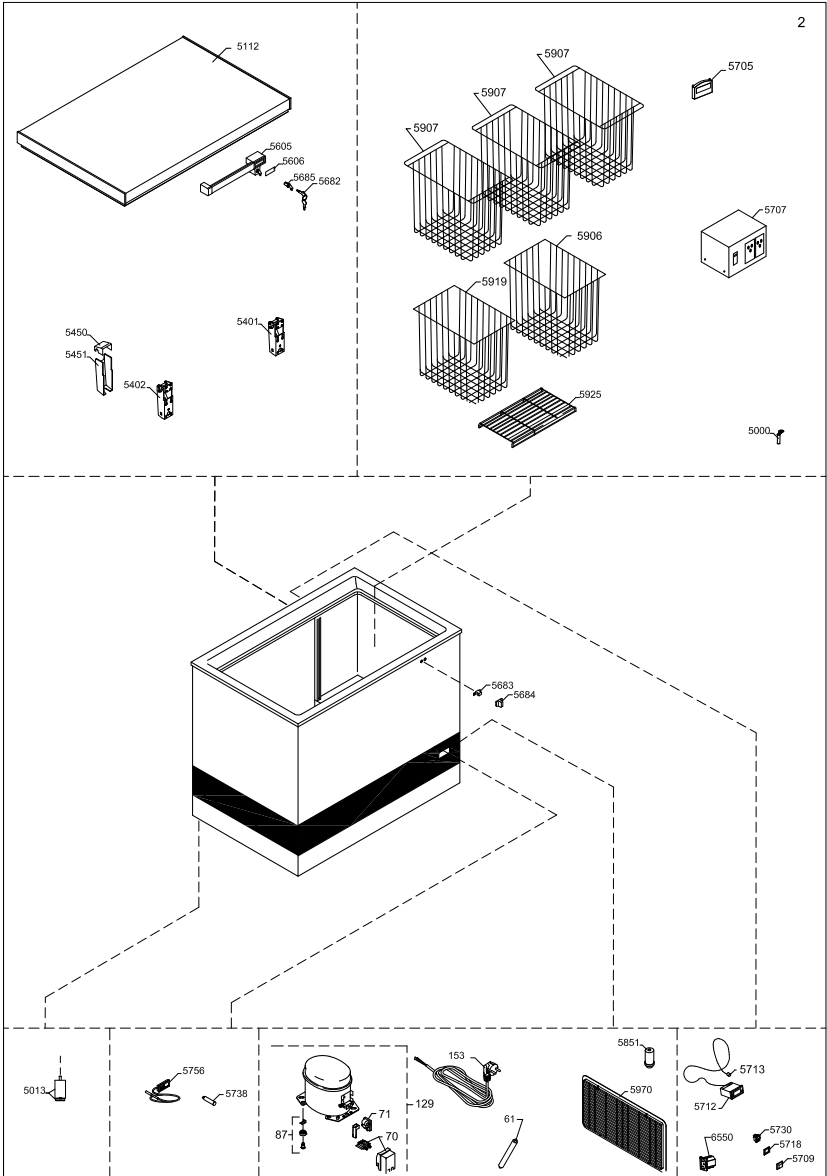
Complete spare part list VLS354A AC

<u>Position</u>	<u>Item number</u>	<u>Item name</u>
0061	0-6538001	Filter drier,
0070	A921001	Cover + wiring clamp
0087	0-6038175	Base plate fittings, complete
0129	8-036038255	Compressor - NLE 9KTK , Complete, R600a
0153	7535254	Power cord EU H05VV-F 3G1 C13 jack 90°
5000	3010049	Drain plug
5013	3040400	Adjustable foot
5116	5000911213	Lid foamed without handle/hinges
5402	1510135	Hinge with spring
5450	3011135-01	Top part for hinge cover
5451	3010032-01	Bottom part for hinge cover
5605	304090501	Handle with lock (push and turn)
5606	8090342-94	Inlay for handle,
5682	1510238	Nikel-plated Key
5683	2040145	Catch for handle
5684	3010265-01	Cover for catch
5685	600098801	Lock with keys,(push and turn)
5705	7020406	Temperature monitoring device Fridge-Tag 2
5707	7020248	Voltage stabilizer
5709	7060104	Frame for cover rocker switch bezel
5712	7095514	Electronic controller ERC 111A
5713	7020960	Sensor NTC S1, Cool
5718	7060105	Rocker Switch Cover
5730	7020245	"on-off" switch
5738	7010138	Pilot lamp, green
5756	7020392-03	Thermometer, solar
5810	A921115	Starting device, 103N0021
5851	6520229	Run capacitor
5874	A921004	Cover + wiring clamp
5907	3510015	Basket 236 mm
5918	3520055	Basket, 190 mm
5919	3520056	Basket 225 mm
5924	3510030	Bottom grating
5970	3010308-01	Motor screen
6550	7080143	Connector

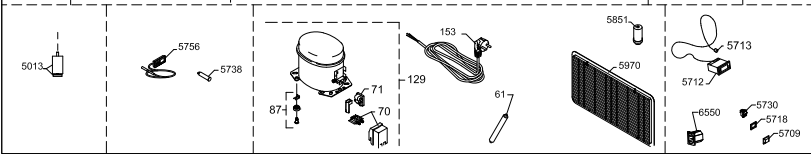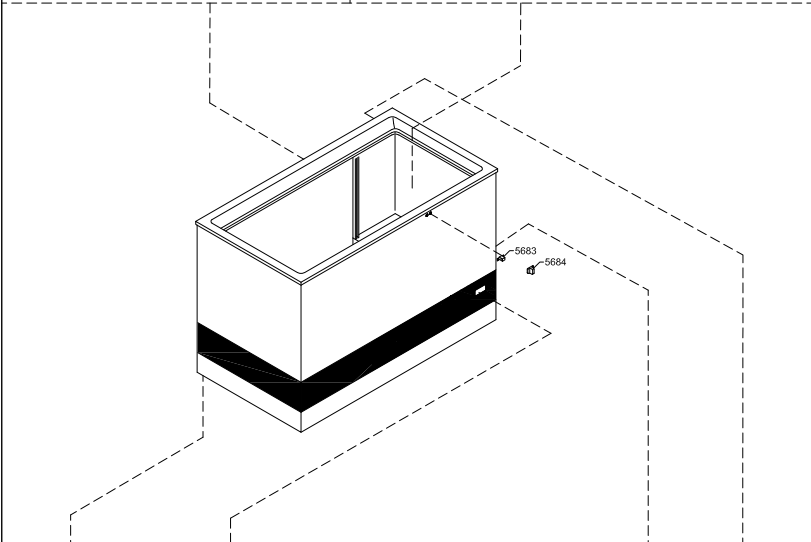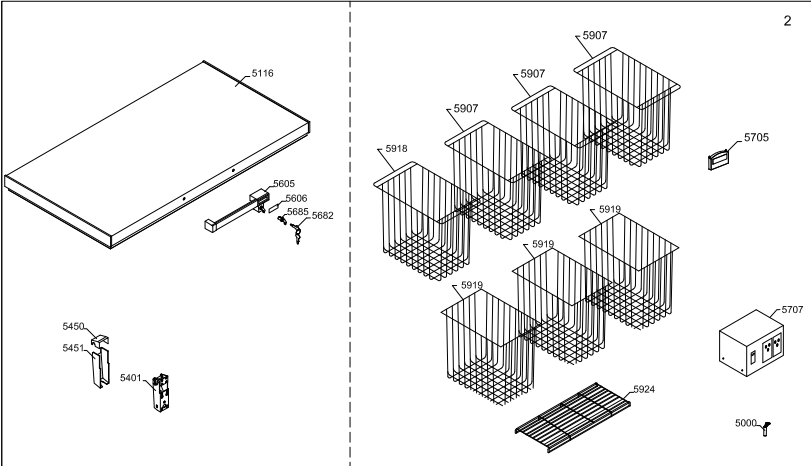Complete spare part list VLS404A AC


Position	Item number	Item name
0061	0-6538001	Filter drier,
0070	A921001	Cover + wirring clamp
0087	0-6038175	Base plate fittings, complete
0129	8-036038255	Compressor - NLE 9KTK , Complete, R600a
0153	7535254	Power cord EU H05VV-F 3G1 C13 jack 90° b
5000	3010049	Drain plug
5013	3040400	Adjustable foot
5118	5000902243	Lid foamed without handle/hinges
5402	1510135	Hinge with spring
5450	3011135-01	Top part for hinge cover
5451	3010032-01	Bottom part for hinge cover
5605	304090501	Handle with lock (push and turn)
5606	8090342-94	Inlay for handle,
5682	1510238	Nikel-plated Key
5683	2040145	Catch for handle
5684	3010265-01	Cover for catch
5685	600098801	Lock with keys,(push and turn)
5705	7020406	Temperature monitoring device Fridge-Tag 2
5707	7020248	Voltage stabilizer
5709	7060104	Frame for cover rocker switch bezel
5712	7095514	Electronic controller ERC 111A
5713	7020960	Sensor NTC S1, Cool
5718	7060105	Rocker Switch Cover
5730	7020245	"on-off" switch
5738	7010138	Pilot lamp, green
5756	7020392-03	Thermometer, solar
5810	A921115	Starting device, 103N0021
5851	6520229	Run capacitor
5874	A921004	Cover + wirring clamp
5906	3510016	Basket 275 mm
5907	3510015	Basket 236 mm
5923	3510028	Bottom grating
5970	3010308-01	Motor screen
6550	7080143	Connector

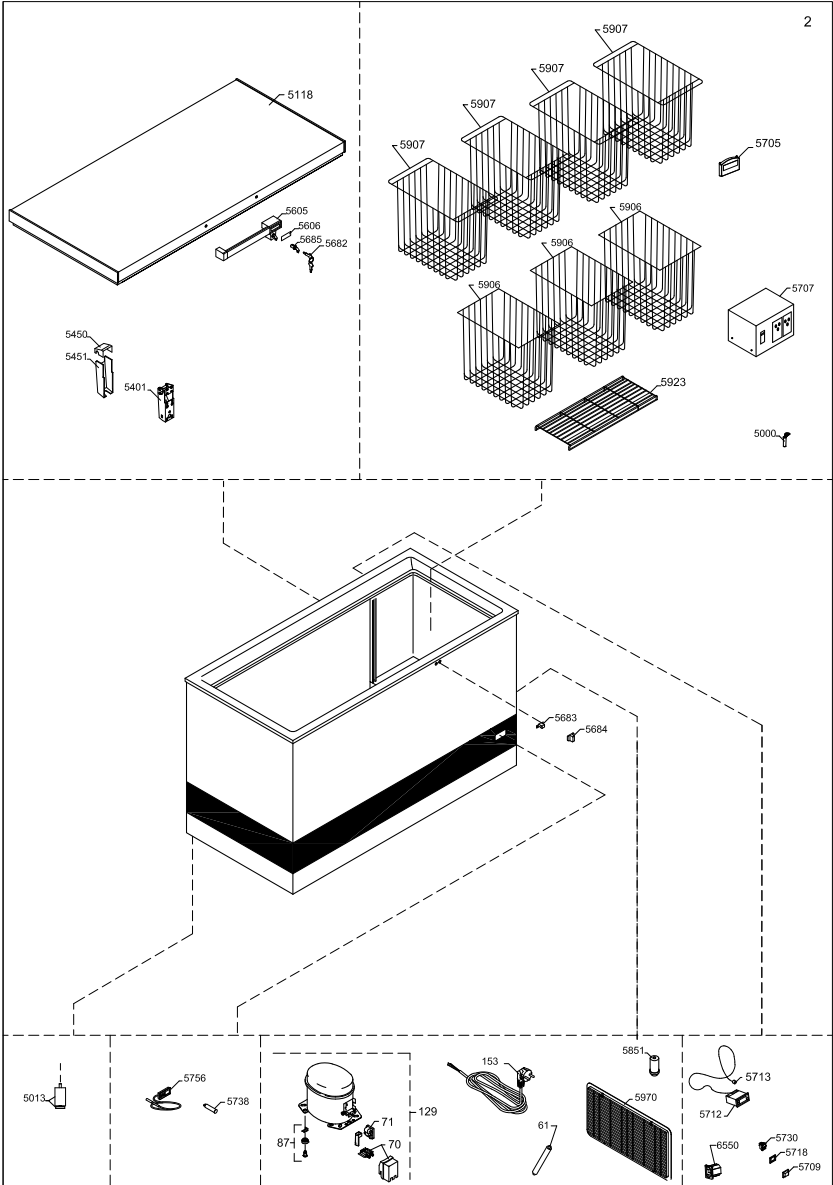Complete spare part list VLS504A AC

<u>Position</u>	<u>Item number</u>	<u>Item name</u>
0061	0-6538001	Filter drier,
0087	0-6038175	Base plate fittings, complete
0129	8-03606510309	Compressor - HXK12AT , Complete,
0153	7535254	Power cord EU H05VV-F 3G1 C13 jack 90°
0550	6520510	Run cap. 4uF/4,8 receptacles
1126	0-A921012	Cover + wirring clamp
1141	6520511	Starting device PTC
2170	0-A9301260103	Distance piece
5000	3010049	Drain plug
5013	3040400	Adjustable foot
5052	2040504	Fittings for condenser
5053	6010068	WOT condenser
5122	5000914293	Lid foamed without handle/hinges
5402	1510136	Hinge with spring
5450	3011135-01	Top part for hinge cover
5451	3010032-01	Bottom part for hinge cover
5605	304090501	Handle with lock (push and turn)
5606	8090342-94	Inlay for handle,
5682	1510238	Nikel-plated Key
5683	2040145	Catch for handle
5684	3010265-01	Cover for catch
5685	600098801	Lock with keys,(push and turn)
5705	7020406	Temperature monitoring device Fridge-Tag :
5707	7020248	Voltage stabilizer
5709	7060104	Frame for cover rocker switch bezel
5712	7095514	Electronic controller ERC 111A
5713	7020960	Sensor NTC S1, Cool
5718	7060105	Rocker Switch Cover
5730	7020245	"on-off" switch
5738	7010138	Pilot lamp, green
5756	7020392-03	Thermometer, solar
5870	7030309	Fan
5906	3510016	Basket 275 mm
5907	3510526	Basket 236 mm
5917	3510525	Basket 117 mm
5931	3510523	Bottom grating
5970	3010308-01	Motor screen
6550	7080143	Connector

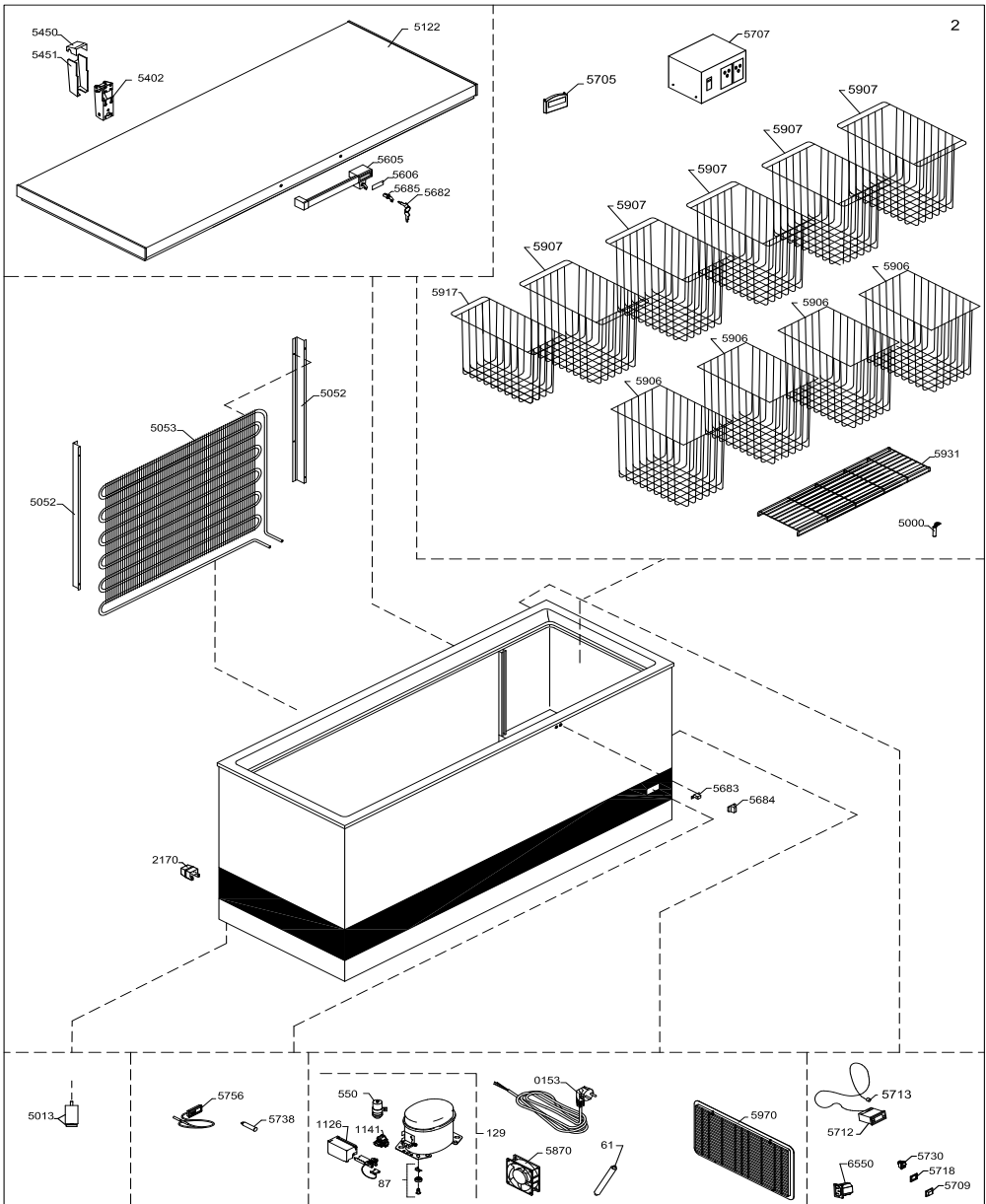
Vital components

Position Item no Description


VLS204AAC & VLS304AAC:
0129 8-036038297 Compressor

VLS354AAC & VLS404AAC:
0129 8-036038255 Compressor

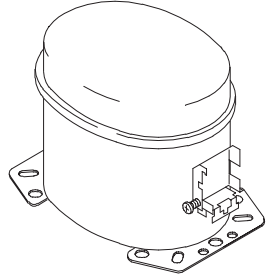VLS504AAC
0129 8-03606510309 Compressor

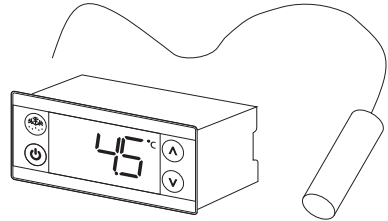
5713 7020961 Sensor
5712 7095514 Thermostat


VLS204,304,354,404AAC
5851 6520229 Run capacitor

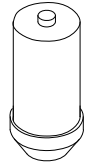VLS504AAC
0550 6520510 Run capacitor

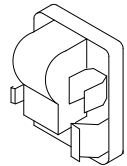
VLS204,304,354,404AAC
0071 A921115 Starting device

VLS504AAC
1141 6520511 Starting device

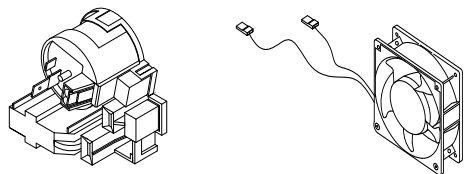5870 7030309 Fan

Health and safety guidance – Warning!

Before any repair job be aware of following!

WARNING:

Before servicing or cleaning the appliance, disconnect it from power source.

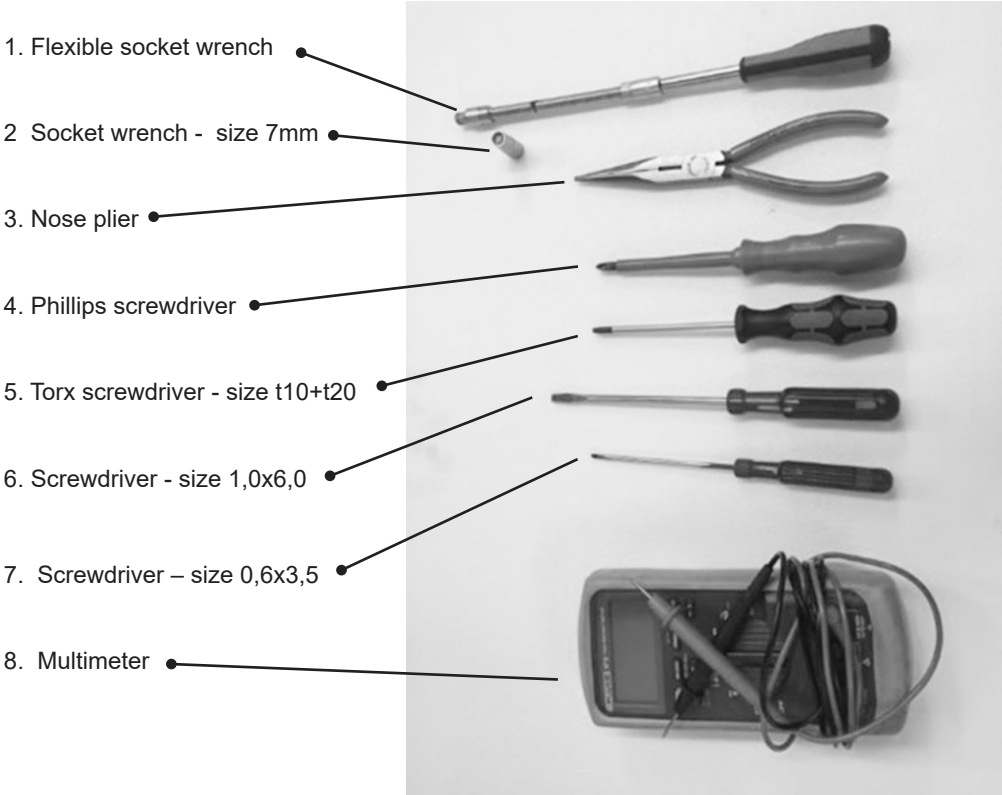
WARNING:

Danger risk of fire or explosion. Flammable refrigerant used. To be repaired only by trained personnel.

Required basic tools


Proposed additional service kit/items

Sealing kit

Tar tape


Extra self-tapping screws

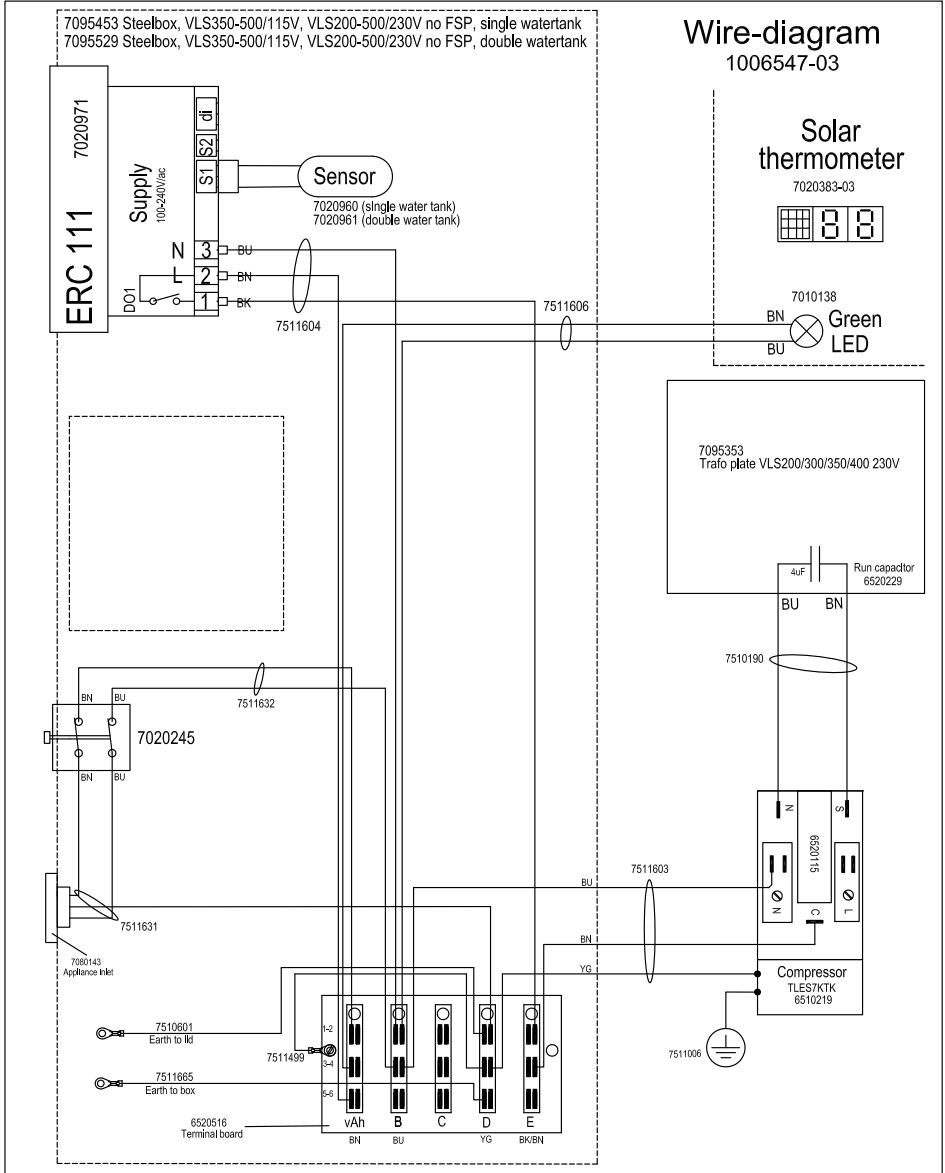Wire diagram VLS204A AC / 304A AC


Wire diagram VLS354A AC / 404A AC


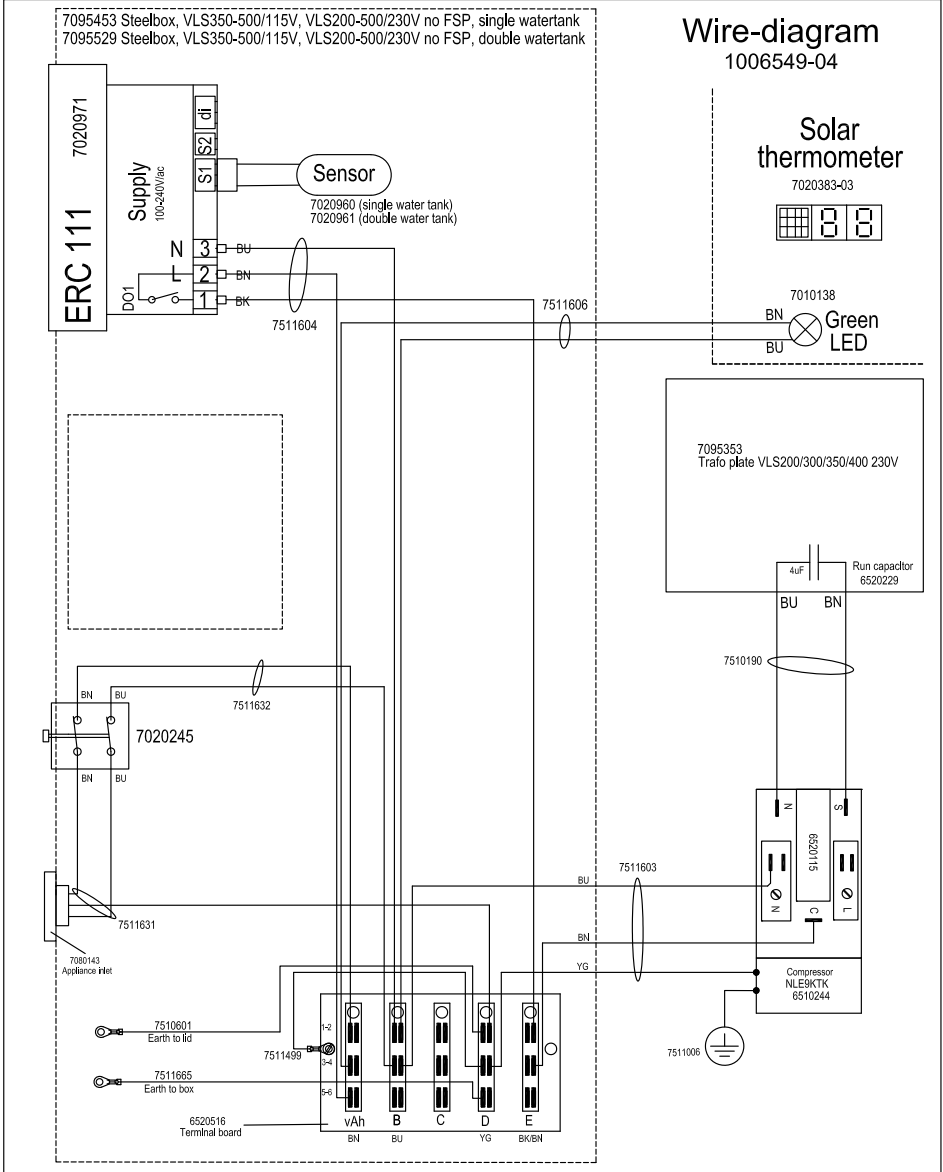
Wire diagram VLS504A AC

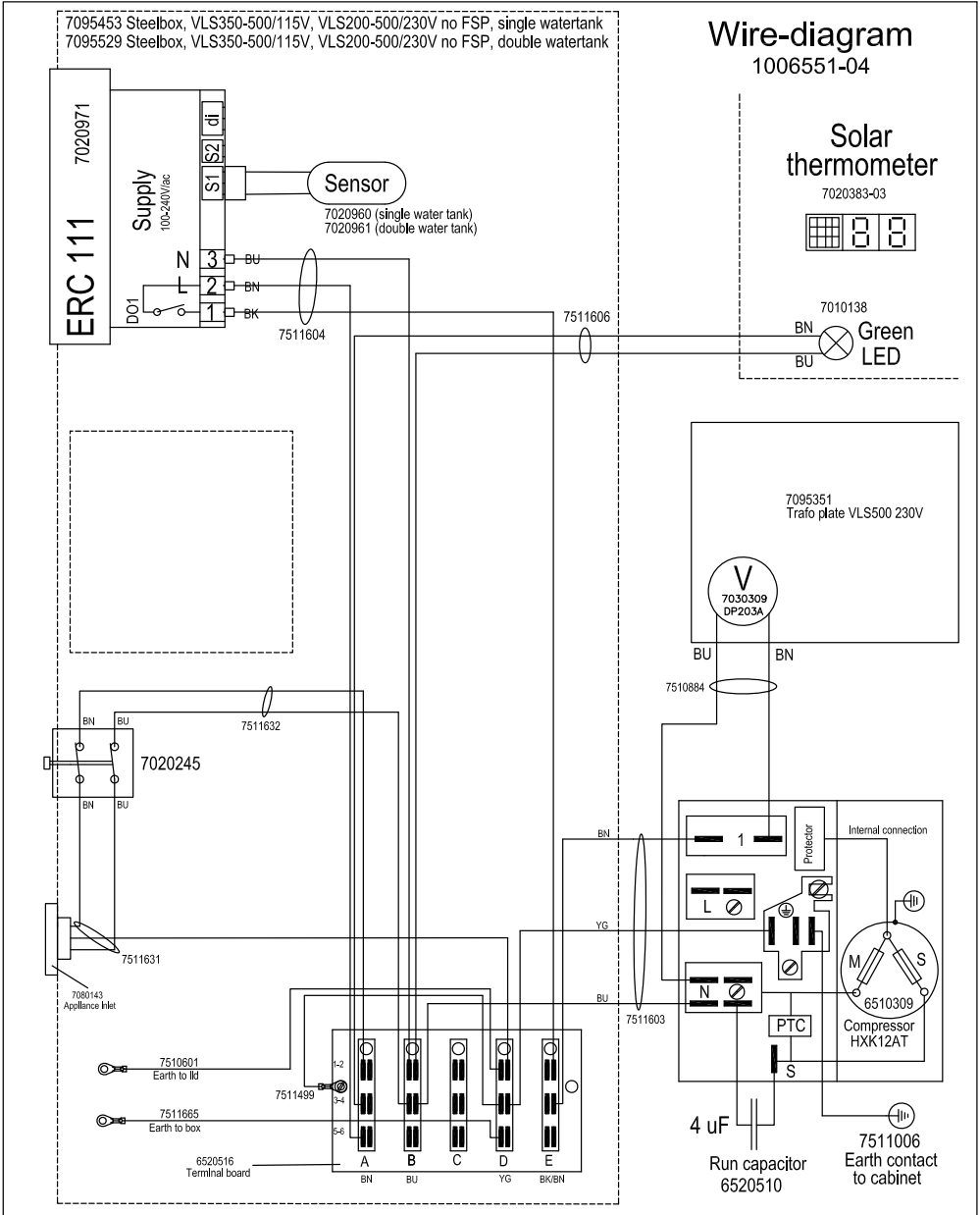
Motor compartment

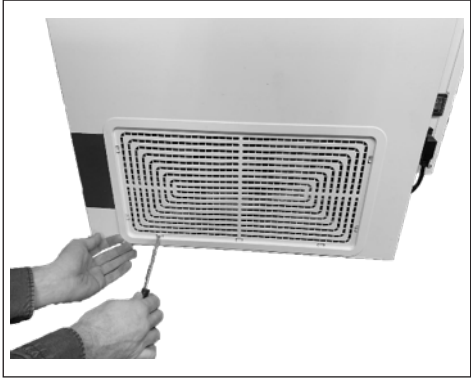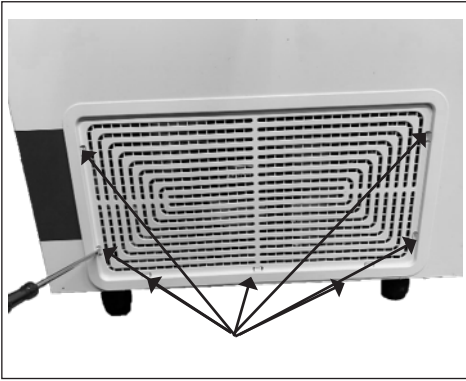How to get access to the motor compartment.

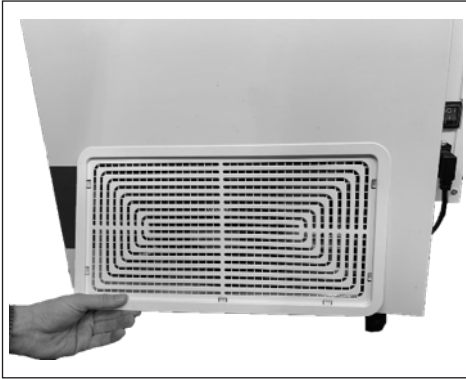1. Use a screwdriver to unlock all 7 clamps.

2. Unlock all 7 clamps

3. Gently pull the compressor grille.

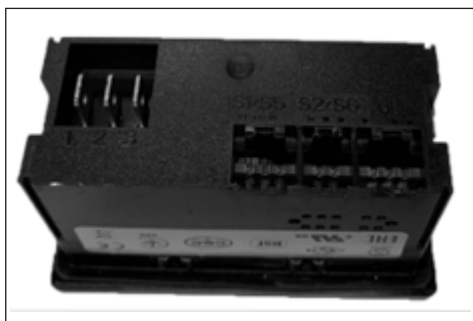
Thermostat replacement

The thermostat is in the junction box located on the rear of the appliance

Frotal con pantallaFront with Display


Side view.


Back with 3x sensor-wire sockets

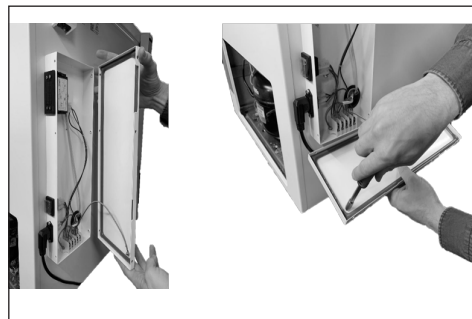1: Open the junction box by unscrewing the 6 Torx 20 screws and unmount lid

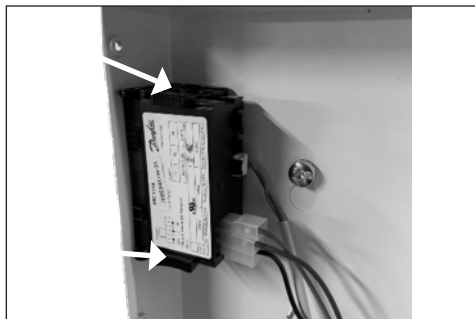
2: Remove grounding cable from junction box by unscrewing bolt with a wrench 6mm

3: Remove 2 clamps from thermostat

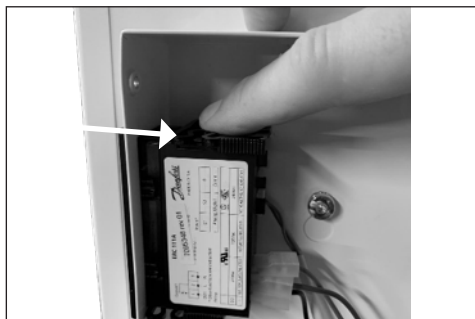Fixing clamp


4: Use your finger to press and slide the lower clamp backwards to remove from the thermostat body


5: Use your finger to press and slide the lower clamp backwards to remove from the thermostat body

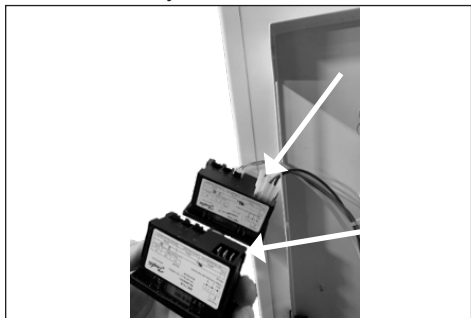6: Push the thermostat out

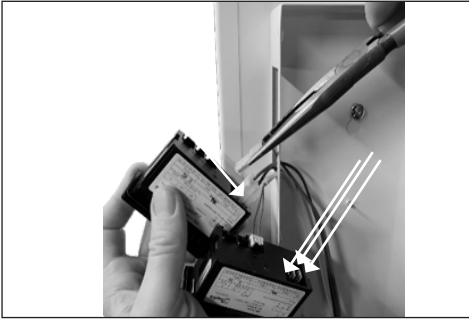7: Switch the sensor wire plug from old socket to the new

8: Switch the 3 power the wires from old to new thermostat one by one

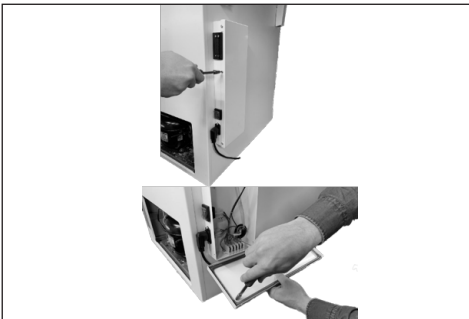
9: Put the thermostat back in place

10: Use your finger to press and slide the upper clamp back in place to secure fixture of the thermostat

11: Use your finger to press and slide the upper clamp back in place to secure fixture of the thermostat

12: Remount grounding wire and junction box lid

Thermostat adjustment

The thermostat is default factory set, and as a general rule not supposed to be adjusted.

In the event, it is assessed by a cold chain technician that a thermostat adjustment is required, follow this instruction.

Steps:

1. Press < > "up/down" and hold 5 seconds to access the menu.
2. Press on/off button  x 2 times till you see numbers flashing
3. Push either < for adjusting lower or > for adjusting higher
4. To save press on/off button  x 1 time then freezing symbol  2 time

Warning!

Thermostat tampering can lead to unsatisfactory cooling, risking damage to stored vaccines.

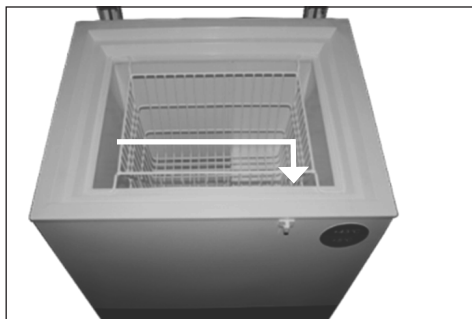
If adjustment is required ONLY to be performed by trained cold chain technicians.

Adjust the controller max. 1°C at a time.


After adjustment monitor appliance carefully for min 24 hours


Thermostat sensor replacement

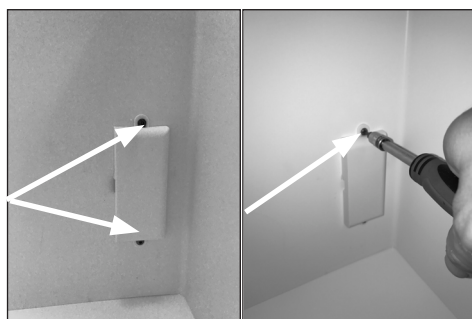
The thermostat sensor is placed inside the compartment of the appliance.

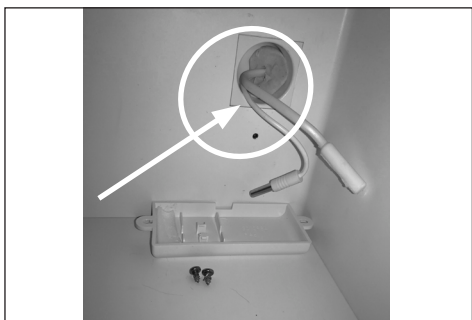
1: Dismount the sensor cover by loosening the 2 x torx screws – size 10.

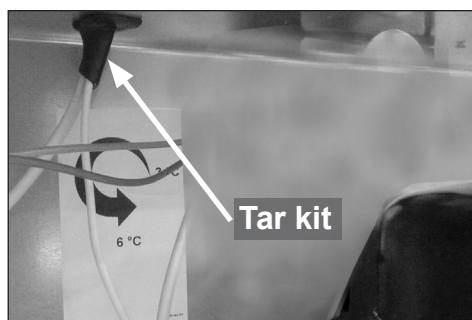
2. Remove the wire sealing and the sensor from the cover.

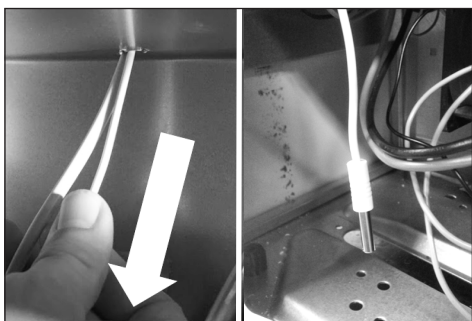
3. Remove black sealing tar-kit and gently pull the white wire until the probe is visible.

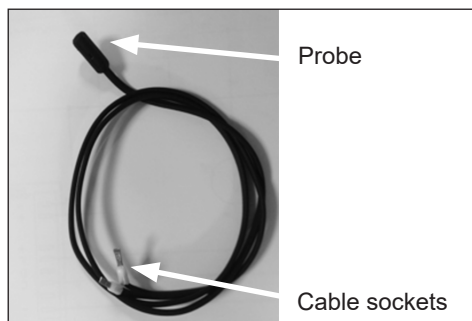
4. Gently pull the white wire until the probe is visible.

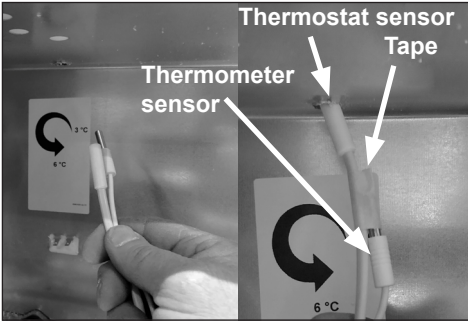
5. The thermostat sensor comes with probe, wire and cable socket.

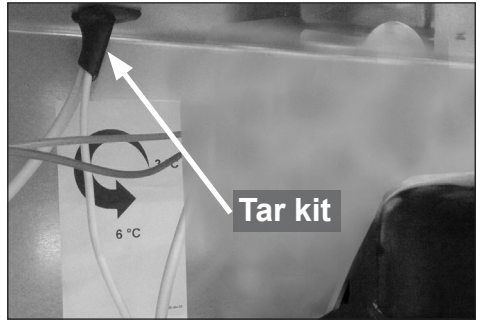6. When re-mounting the new thermostat sensor, use removable tape.

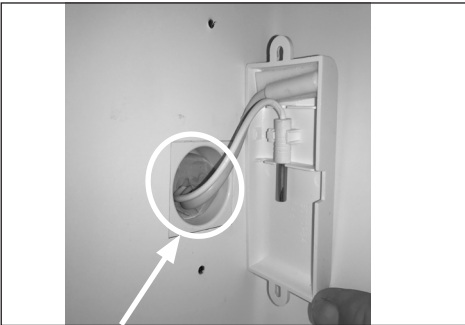7. **IMPORTANT!** When re-mounting the new thermostat sensor remember to properly seal the wire feed through.

8. **IMPORTANT!** When re-mounting the new thermometer make sure the wire sealing plug is placed properly.

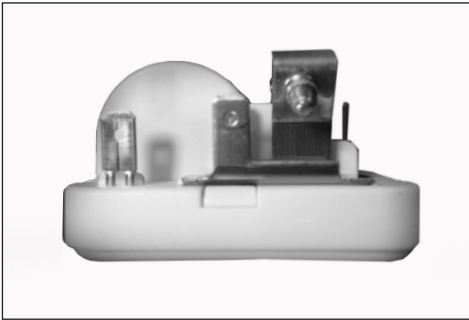
Starting device replacement for VLS204 to 404A AC

Front with terminals.

Back with connection plug.


Side view.

The starting device is mounted on the left side of the compressor.

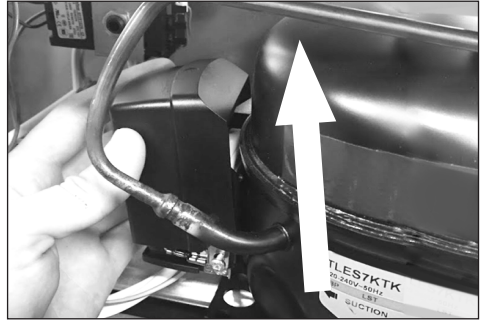
1. Dismount the cover for starting device by loosen the phillips screw.


2. Push the plastic cover up.

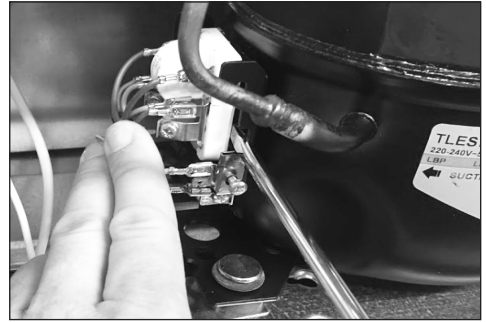
3. Pull the cover back to loosen.

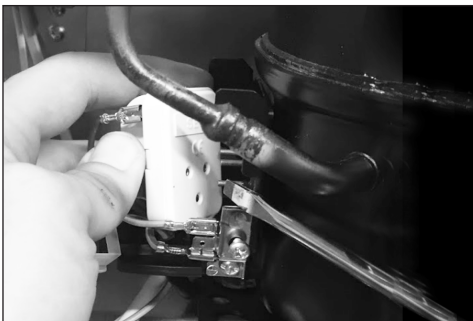
4. Use a screw driver and gently remove the starter from the socket of compressor.

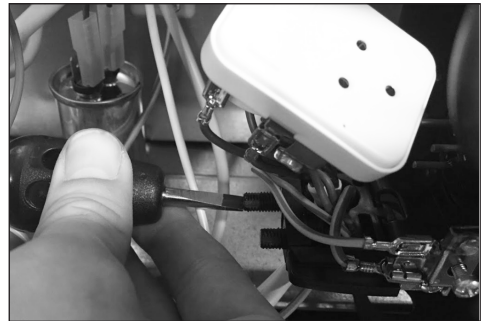5. Starting device loose from socket

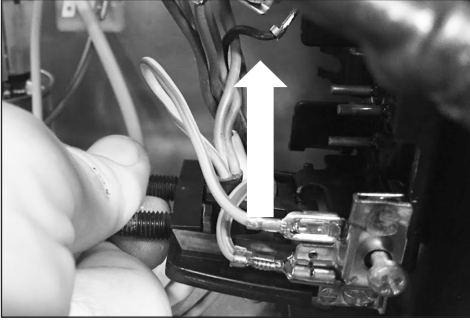
6. Use a small screwdriver or a wrench size 6 – to unmount the wire fastener

7. Push the plastic bracket up.

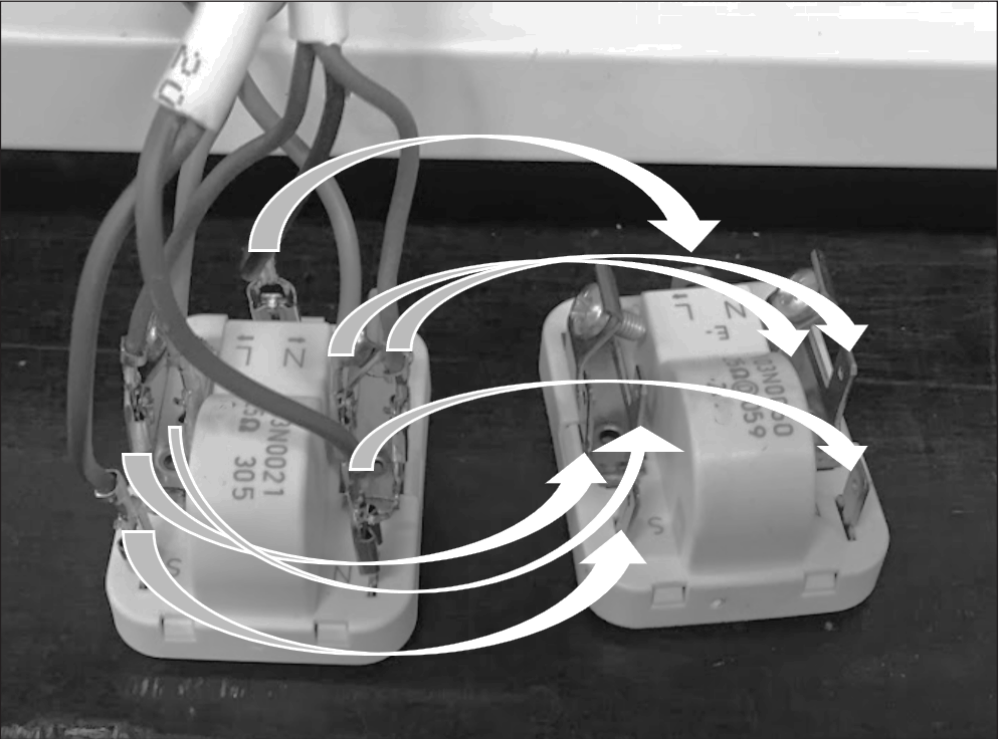
8. Use a nose plier to unmount the wire sockets from starting device.


9. Exchange the wires 1/1 from the old starting device to the new one.

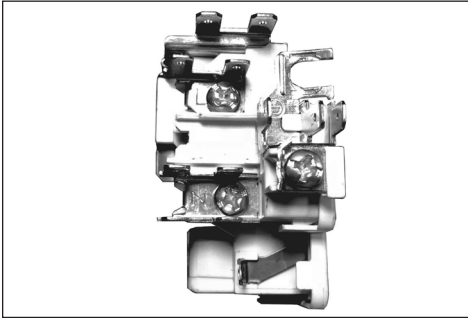
Starting device replacement for VLS504A AC

Front with terminals.

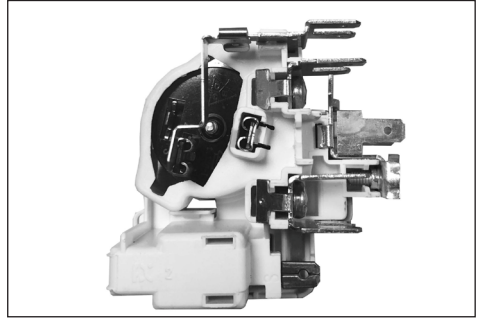
Right side view

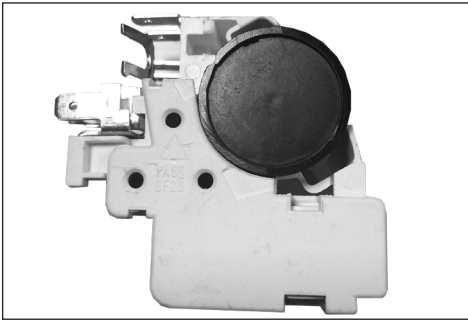Left side view.

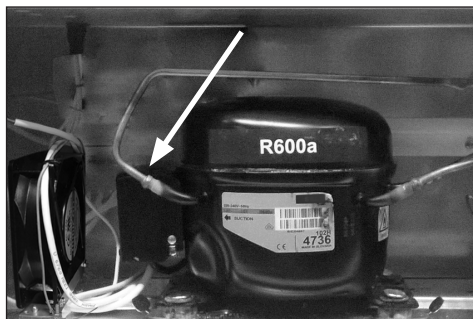The starting device is mounted on the left side of the compressor.


The starting device is placed in the left side of the compressor compartment, fixed to the compressor bracket.

1. Dismount the cover by pushing the cover lock from right to the left with a flat screw driver.

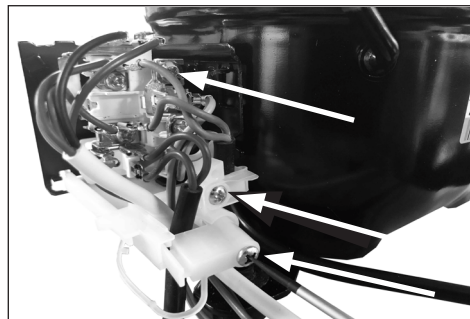
2. Pull out the cover to remove


3. Use a screw driver to loosen 3 x screws.

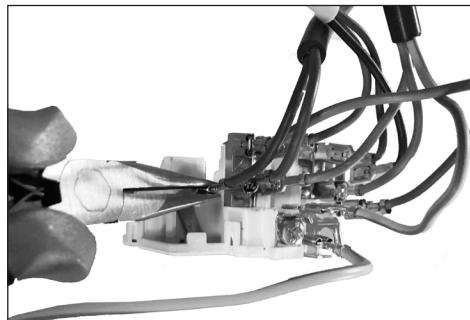4. Use a screwdriver to loose the starting device from socket and pull.

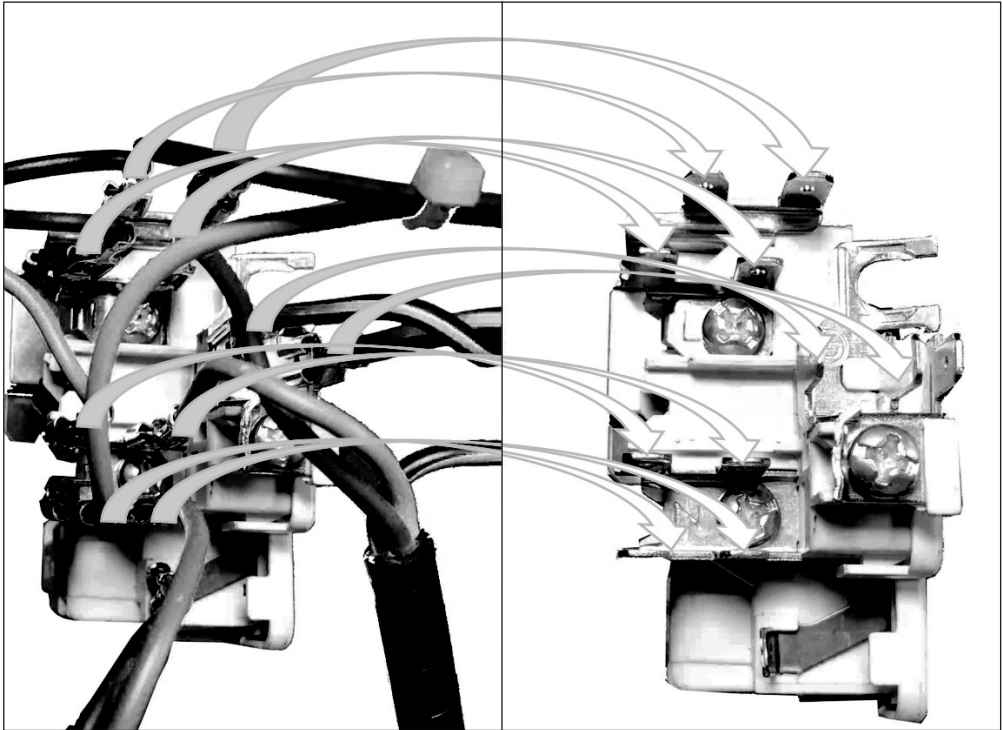5. Use a nose plier to unmount the wire sockets from starting device.

6. Exchange the wires 1/1 from the old starting device to the new one.

Run capacitor replacement VLS204 to 404A AC

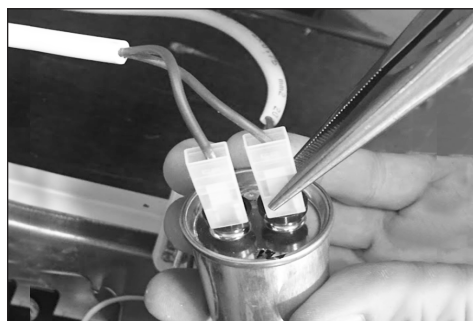
The run capacitor is placed in the left side of the compressor compartment, fixed to the terminal bracket.

1. Remove nuts to unmount assembly plate for run capacitor. Lift up and pull out the assembly plate.

2. Loosen the nut from the assembly plate by using a wrench M13.

3. Unmount the 2 wires by using a nose plier.

4. Exchange the wires 1/1 from the old capacitor to the new one

5. When remounting assembly plate remember remounting of earth wires.

Run capacitor replacement for VLS504A AC

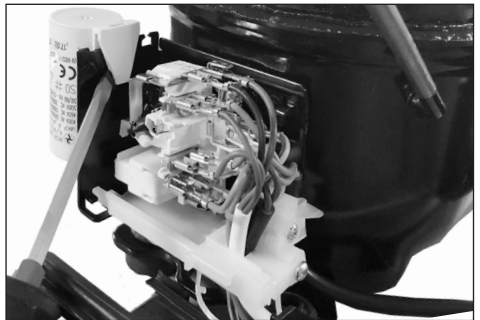
The run capacitor is placed in the left side of the compressor compartment, fixed to the compressor bracket.


1. Dismount the cover by pushing the cover lock from right to the left with a flat screw driver.


2. Pull out the cover to remove

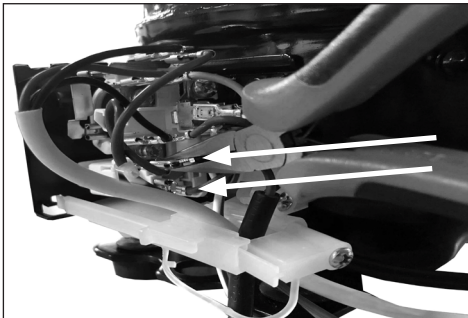3. Flip out the clamp to free the capacitor from the bracket.

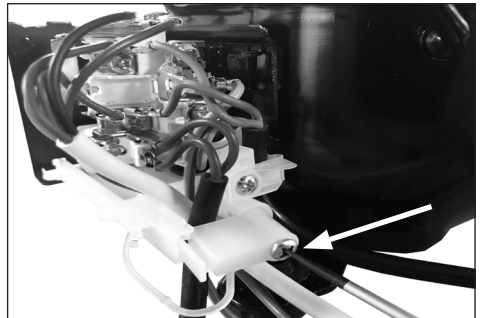4. Use a nose plier to unmount the wire sockets.


5. Loosen the screw for cord relieve and pulle out the wire. The capacitor can be remounted in reverse order.


Thermometer replacement

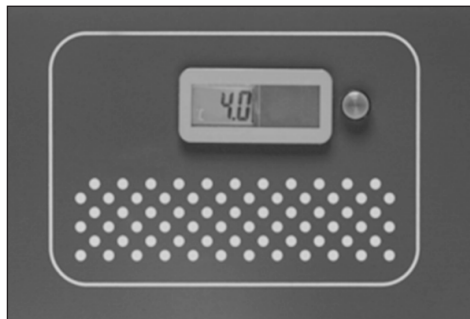
The thermometer is placed in at the front of the appliance.


Thermometer display.

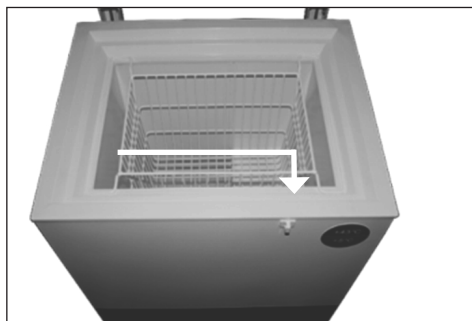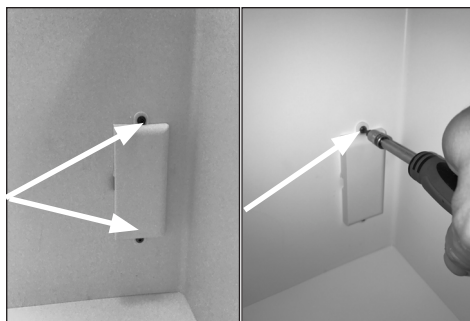The thermometer sensor is placed inside the compartment of the appliance.

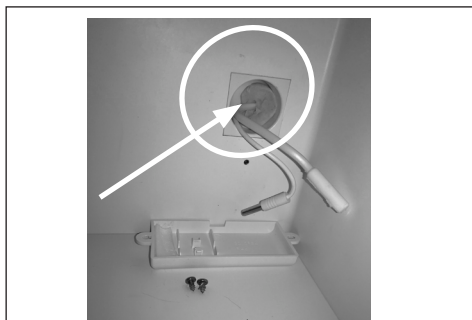
1: Dismount the temperature sensor cover by loosen the 2 x torx screws – size 10.

2. Remove the wire sealing and the sensor from the cover.

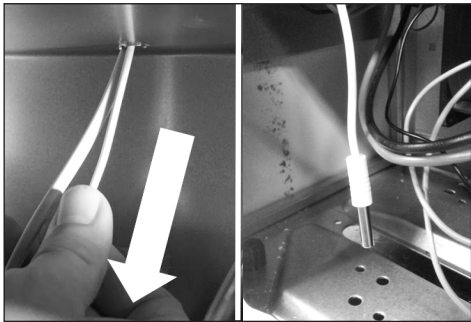3. Remove black sealing tar-kit and gently pull the white wire until the probe is visible.


4. Gently pull the white wire until the probe is visible.

5. Use a screw driver to gently push the socket of the thermometer.

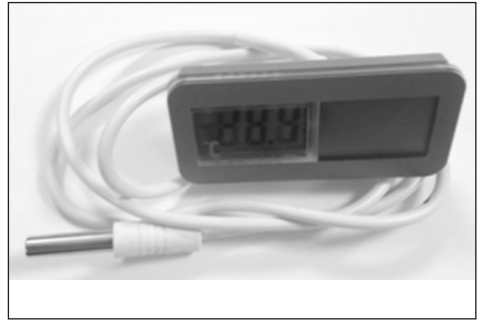
6. Temperature monitor is loose from cabinet.


7. Thermometer comes with wire and sensor.

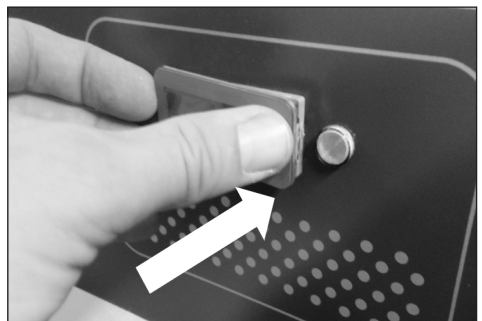
8. Installation of thermometer display.

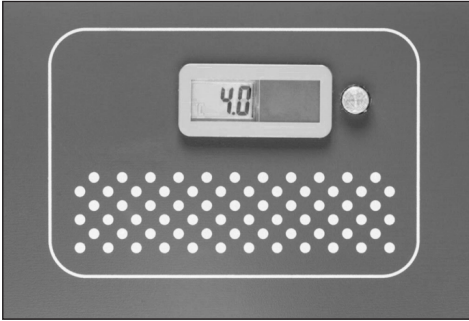9. Push until display is fixed to cabinet.

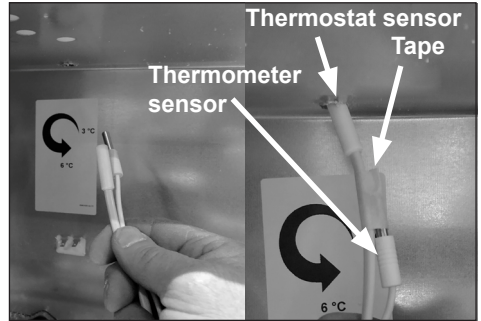
10. Thermometer display is in place.

11. When re-mounting the new thermometer sensor, use removable tape.

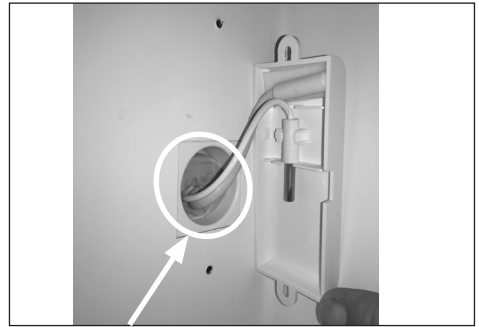12. **IMPORTANT!** When re-mounting the new thermometer remember to properly seal the wire feed through.


13. **IMPORTANT!** When re-mounting the new thermometer make sure the wire sealing plug is placed properly.


Fan replacement VLS504A AC

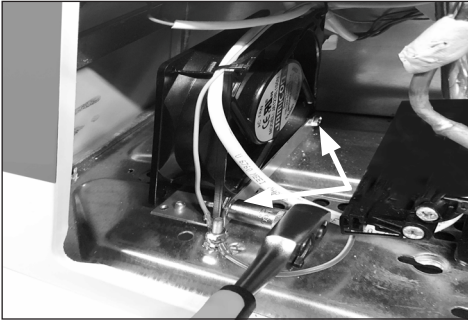
The fan is placed at the left side of the compressor.

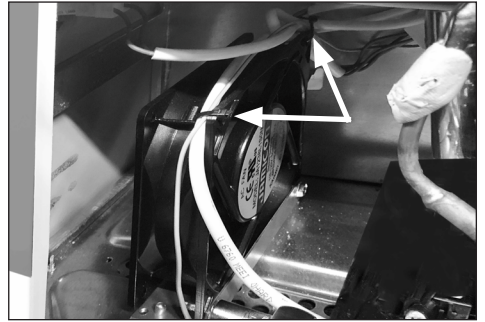1. Dismount the fan from the bracket by losen the 2 hexagon bolts with a Socket wrench.

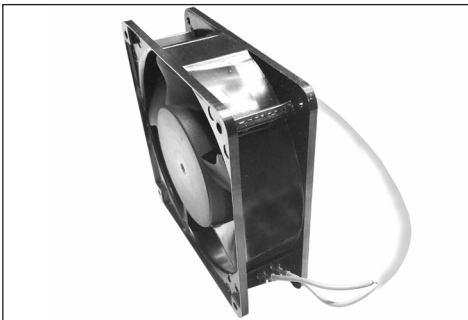2. Remove strips for free access to the fan. When mounting the fan please remember to secure the wires to the fan again.

3. Use the nose plier to unmount the 2x cable sockets.

Compressor replacement

Procedure of compressor switch.

1: **WARNING!** Drain coolant R600a from refrigeration system by vacuum suction.

2: **IMPORTANT!** Blow refrigeration system with NO/Nitrogen

3: Cut

A: Suction and pressure tube

B: Capillary tube

C: Dry filter

4: Dismount starting device

5: Dismount old compressor

6: Insert new compressor

7: Solder

A. Suction and pressure tube

B. Capillary tube

C. Dry filter

8: Install starting device

IMPORTANT! When solder copper tubes to iron tubes use silver tin

Filling of new refrigerant

9: Drain refrigeration system by vacuum suction

10: Check type sticker for required amount of R600a to fill on refrigerant system

On-site checklist

- Is the green diode in the control panel on (Power check)?
- Temperature records (manual records, FT2 data)
- Is the internal temperature inside the acceptable range of +2° to +8°?
- Is the vaccine compartment clean and without condensation (water)?
- Is the Compressor running?
- Are baskets used and in place?
- Is the appliance placed according to instructions in the manual?
- Does the lid close tight to the cabinet and is the lid gasket in good condition?
- Is the grille for the compressor compartment clean?
- Are all electrical components working properly?
- Overall condition of the cabinet – internal and external: any corrosion, rusting, cracks?
- Inspection of the refrigeration system (the condenser, evaporator, the whole refrigeration circuit/line)?


Trouble shooting

Fault	Possible cause	Remedy
Compressor is not running.	<p>Be patient, it is most likely that the compressor will start within a few minutes.</p> <p>Power for appliance is not plugged in</p> <p>Power for stabilizer is not plugged in</p> <p>Power for appliance is not switched on</p> <p>Power for stabilizer is not switched on</p> <p>Stabilizer power fuse has blown</p> <p>Stabilizer 6 min. time delay is in progress</p>	<p>If this is not the case, check the following:</p> <p>Check that the power for appliance is plugged in</p> <p>Check that the power for stabilizer is plugged in</p> <p>Check that the power for appliance is switched on</p> <p>Check that the input voltage is within the input range of stabilizer</p> <p>Check fuse</p> <p>Wait for the delay to end</p> <p>- If the above is OK, call technical supervisor.</p>
Compressor is running, and the temperature is too high.	<p>The ventilation grille is blocked.</p> <p>The lid is not closed properly.</p> <p>The temperature in the room in which the appliance is installed is too high.</p>	<p>Ensure unhindered air circulation.</p> <p>Ensure that the lid is closed properly.</p> <p>Shield the appliance against direct sun light and ensure more ventilation to the room.</p>
No temperature is displayed.	<p>The Thermometer is broken.</p> <p>There is not enough light for the solar sensor.</p>	<p>Change the thermometer.</p> <p>Turn on the light.</p>

Technical support

When contacting Vestfrost Solutions technical support please supply below information:

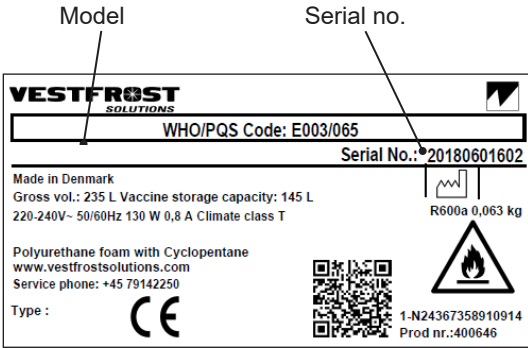1. Model
2. Serial number
3. What is the issue

Contact:

Vestfrost Solutions

Tel. +45 75142250

cce-service@vestfrostolutions.com

Or visit our service-center webpage:

<http://www.vestfrostolutions.com/service-center/>

Recycling procedures

Information for Users on Collection and Disposal Old Equipment and used Batteries

This symbol on the products, packaging, and/or accompanying documents mean that used electrical and electronic products and batteries should not be mixed with general household waste. For proper treatment, recovery and recycling of old products and used batteries, please take them to applicable collection points, in accordance with your national legislation and the Directives 2012/19/EU and 2006/66/EC.

By disposing of these products and batteries correctly, you will help to save valuable resources and prevent any potential negative effects on human health and the environment which could otherwise arise from inappropriate waste handling.

For more information about collection and recycling of old products and batteries, please contact your local municipality, your waste disposal service or the point of sale where you purchased the items.

Penalties may be applicable for incorrect disposal of this waste, in accordance with national legislation.

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