Cabinet WFG48 - WFG49

GB Instructions for use
DK Brugsanvisning
NO Bruksanvisning
SE Brugsanvisning

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Warning

WARNING:

Keep ventilation openings in the appliance's cabinet or in the built-in structure clear of obstruction.

WARNING:

Do not use other **mechanical devices** or other means to accelerate the defrosting process than those recommended by the manufacturer.

WARNING:

Do not damage the refrigerant circuit.

WARNING:

Do not use **electrical appliances** inside the storage compartment, unless they are of a type recommended by the manufacturer.

WARNING:

Danger: Risk of fire or explosion if flammable refrigerant is used. To be repaired only by trained personnel.

WARNING:

When positioning the appliance, ensure the power cord is not trapped or damaged.

WARNING:

Do not locate multiple portable socket-outlets or portable power supplies at the rear of the appliance.

- Always keep the keys in a separate place and out of reach of children.
- Children should be supervised to ensure that they do not play with the appliance.
- This appliance is not intended for use by persons (including children) with reduced physical, sensory or mental capabilities, or lack of experience and knowledge, unless they have been given supervision or instruction concerning use of the appliance by a person responsible for their safety.
- Before servicing or cleaning the appliance, unplug the appliance from the mains or disconnect the electrical power supply.
- If the supply cord is damaged, it must be replaced by the manufacturer, its service agent, or similarly qualified persons in order to avoid a hazard.

Contents

Warning	2
Before use	4
Technical data.....	4
Get to know your cooler.....	5
Installation and start-up	6
Savings in energy consumption in the use ..	6
Operation and function	7
Change of hinge side	9
Defrosting, cleaning, and maintenance	12
Fault finding	13
Warranty, spare parts and service	14
Disposal.....	15
CE Declaration of conformity.....	Page A

- Frost formation on the interior evaporator wall and upper parts is a natural phenomenon. Therefore, the appliance should be defrosted during normal cleaning or maintenance.
- Please note that changes to the appliance construction will cancel all warranty and product liability.
- This appliance is intended to be used exclusively for the storage of bottles or cans.
- This appliance is intended to be used in household and similar applications.
- Do not store explosives, such as aerosol cans with flammable propellants in the unit.

CLASS 1 LED PRODUCT

The instructions shall include the substance of the following:

This appliance is intended to be used in household and similar applications such as

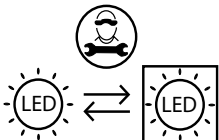
- staff kitchen areas in shops, offices and other working environments;
- farm houses and by clients in hotels, motels and other residential type environments;
- bed and breakfast type environments;
- catering and similar retail and non-retail applications.

Before use

Before operating your new appliance, please read the following instructions carefully as they contain important information on safety, installation, operation and maintenance. Keep the instructions for future reference. On receipt, check to ensure that the appliance has not been damaged during transport. Transport damage should be reported to the local distributor before the beverage cooler is put into use.

Remove the packaging. Clean the inside of the cabinet using warm water with a mild detergent. Rinse with clean water and dry thoroughly (see cleaning instructions). Use a soft cloth. If the appliance has been laid down during transport, or if it has been stored in cold surroundings (colder than +5°C), it must stabilise in an upright position for at least one hour before being switched on.

Replaceable light source by a professional

Replaceable control gear by a professional

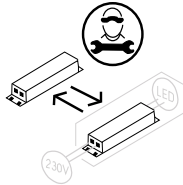
fig. 1

Technical data

Wiring and connections in power supply systems must be all applicable (local and national) electrical codes. Consult these codes lengths and sizes prior to cabinet installation.

This device complies with relevant EU directives including Low Voltage Directive 2014/35/EU and Electromagnetic Compatibility Directive 2014/30/EU

The socket should be freely accessible.

Connect the appliance only to 220/240 V / 50Hz alternating current via a correctly installed earthed socket.

The socket must be fused with a 10-13 A fuse.

If the appliance is to be operated in a non-European country, check on the rating plate whether the indicated voltage and current type correspond to the values of your mains supply.

Information regarding voltage, current or power are given on the rating plate

The power cord may be replaced by a technician only.

The rating plate provides various technical information as well as type and serial number.

Get to know your cooler

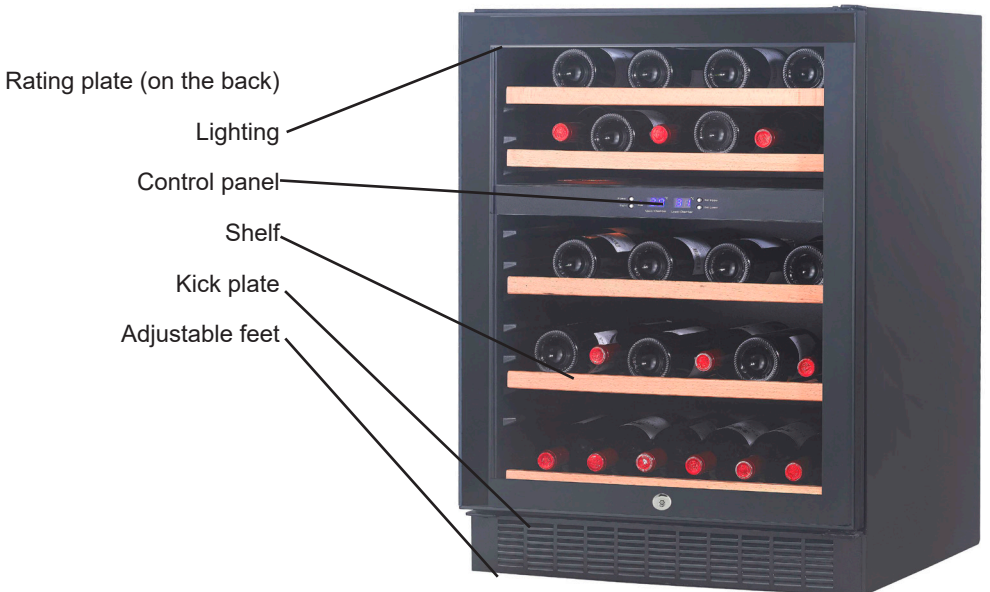
fig. 2

Installation and start-up

Placement

For safety and operational reasons, the appliance must not be installed outdoors.

The appliance should be placed on a level surface in a dry, well ventilated room (max. 75% relative air humidity). Never place the appliance close to sources of heat such as cookers or radiators, and avoid placing it in direct sunlight.

Installation

The surface on which the appliance is to be placed must be level. Do not use a frame or similar. It is important that the appliance be absolutely level. It can be levelled by screwing the adjustable feet at the front of the appliance up or down (fig. 2).

Use a spirit level to check that the appliance is absolutely level sideways.

Along with your new cabinet is a replaceable kick plate which can be fixed with the two screws. These two screws are located in the accessories. Hinge side requires a 5 mm space.

NOTE: The kick plate should not be installed if the unit installed behind a kitchen shelf.

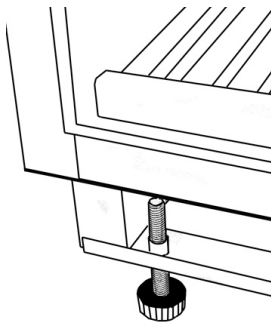
fig. 3

Room temperature

The climate class is stated on the rating plate (see page 11).

Note: The appliance operates best at room temperatures from 10°C to 32°C.

The climate class is shared on the name-plate. This specifies the optimum room temperature.

Climate class	Optimum room temperature
SN	+10 °C to +32 °C
N	+16 °C to +32 °C
ST	+16 °C to +38 °C
T	+16 °C to +43 °C

Max. 19 kg

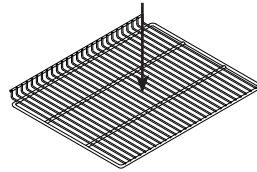
fig. 4

Savings in energy consumption in the use

The energy consumption reduces ..

- 1)..the lower the ambient temperature is.
- 2)..the warmer thermostat set point is.
- 3)..the lower the number and the shorter the length of the door openings is.

Operation and function

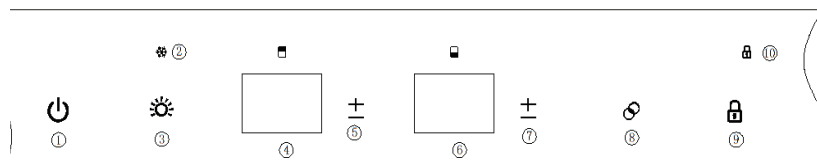
fig. 5

1. **Power ON/OFF button (beep sound)**
Press this button to turn on the appliance. Press and hold this button for 3 seconds. The left display shows a countdown from 3 to 1, then the appliance turns off.
2. **Indicator light**
Displays compressor in operation.
3. **Light mode button (beep sound)**
There are three light mode types (L1/L2/L3). The default setting is L1 mode.
L1: Door open: light on. Door closed: light off.
L2: The light will stay constantly illuminated whether the door is open or closed. Door open with 100% light brightness. Door closed with 75% light brightness.
L3: Door open: light on. Door closed: light will tune off automatically after 10 minutes.
4. **Left display**
Shows the set temperature of the upper chamber.
5. **Button to adjust upper zone temperature (beep sound)**
+ Each pressure on this button decreases the temperature by one-degree increments. When the setting reaches the lower limit value of 5°C, it will revert to the upper limit value of 10°C. After selecting your desired set temperature, the set value will flash for 5 seconds, after which the display will show the set temperature.
-
6. **Right display**
Shows the set temperature of the lower chamber.
7. **Button to adjust lower zone temperature (beep sound)**
+ Each pressure on this button decreases the temperature by one-degree increments. When the setting reaches the lower limit value of 10°C, it will revert to the upper limit value of 18°C. After selecting your desired set temperature, the set value will flash for 5 seconds, after which the display will show the set temperature.
-

8.  **Light colour button (beep sound)**
Press this button to alternate the light colour. The default setting is white light. Each pressure alternates the light colour between White - Blue - Orange. The button background colour also changes at the same time.
9.  **Child Lock button (beep sound)**
The default setting is unlocked. Press and hold this button for 3 seconds to lock or unlock the appliance. The indicator lamp is on when in locked condition.
10.  **Indicator lamp**
The indicator lamp is on when the appliance is in locked condition.

NOTE

To switch between °C and °F, press and hold the 5 or 7 temperature adjust button for 3 seconds. The display will show the desired setting.

Temperature Alarm

If the inside temperature is higher than 23 °C, the display will show "HI", and after one hour the alarm / buzzer will sound. This indicates that the inside temperature is too high. If the inside temperature is lower than 0 °C, the display will show "LO", and at the same time the alarm malfunction indicator is also active (beep sound).

Open door Alarm

When you forget to close the door, or the door is not closed completely, the alarm / buzzer sounds after 5 minutes. Please close the door completely to avoid loss of the cold air.

Temperature memory function

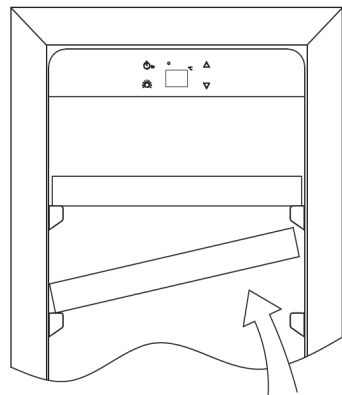
In the event of a power failure, the wine cooler can remember the previous temperature settings, and when the power is recovered, the temperature will revert to the same temperature setting as before the power failure.

Shelves

In order, not to damage the gasket on the door, it is very important to open the door completely before pulling out the shelves.

Shelves should be tilted slightly before being removed or inserted, as shown at the right.

For easy access to the stored content, pull the shelves approximately 1/3 out of the compartment. The wooden shelves are designed with concave grooves, so that tall wine bottles can be stored crossways.

Change of hinge side

The door can be changed from right-hinged to left-hinged and vice versa as follows:

fig. 7.1

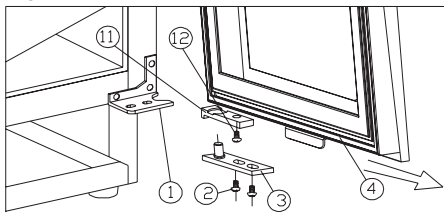
fig. 7.2

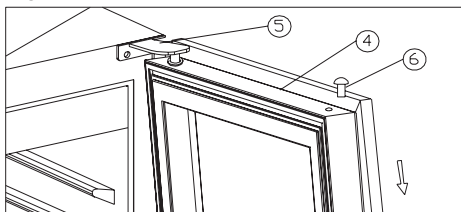
fig. 7.3

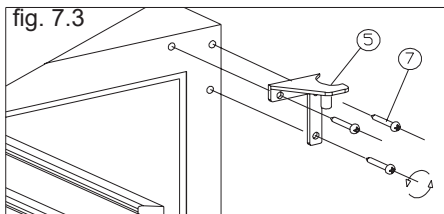
fig. 7.4


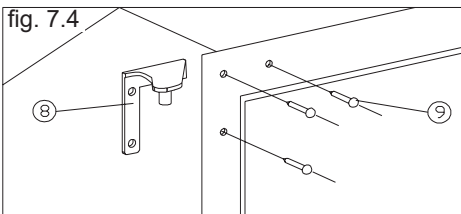
fig. 7.5

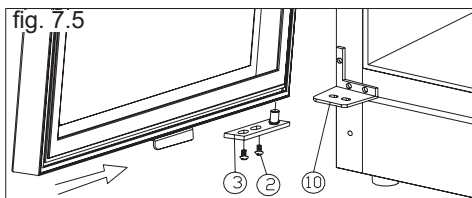Ensure that the cabinet is not connected to the mains and place the cabinet on the back. SUPPORT THE DOOR !

- | | | | | |
|------------------------|------------------|---------------------|---------------|----------------------|
| 1. Lower hinge (right) | 2. Screws | 3. Door hinge | 4. Door | 5. Top hinge (right) |
| 6. Plastic cover | 7. Hinge screws | 8. Top hinge (left) | 9. Hole cover | |
| 10. Lower hinge (left) | 11. Door fitting | 12. Screws | | |

1. Open the door to a 90 degrees angle, take out the 2 screws (2) from lower door hing (3), remove the door (4), Remove the door hinge (3) from the door, (Fig. 7.1).
2. Unscrew the screws (12) from the door fittings (11), and relocate it at the opposite side of the door (fig. 7.1)
3. Remove plastic cover (6) (fig. 7.2).
4. Unscrew the 3 hinge screws (7) from the top hinge (5) and remove it (fig. 7.3).
5. Remove hole covers (9) and mount the spare top hinge (8) at the opposite side of the cabinet with the 3 screws (fig. 7.4).
6. Mount the spare lower hinge (10) (fig. 7.5).
7. Mount and align the door.
8. Insert the hole covers to cover the holes (9).

Automatic door opening function

After the door is closed, press 2-3 seconds on the lower left of the door, after which the push rod starts and sticks out, and the door opens automatically, and then the push rod automatically retracts. After two seconds, the door can be closed manually.

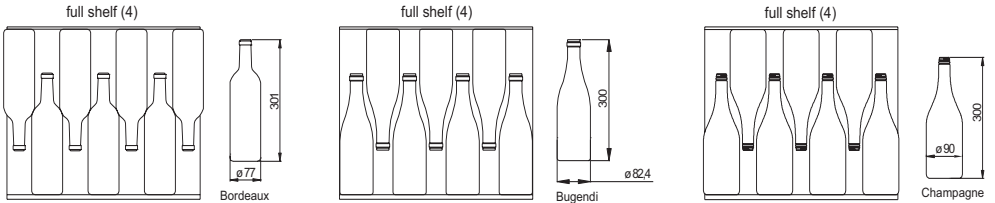
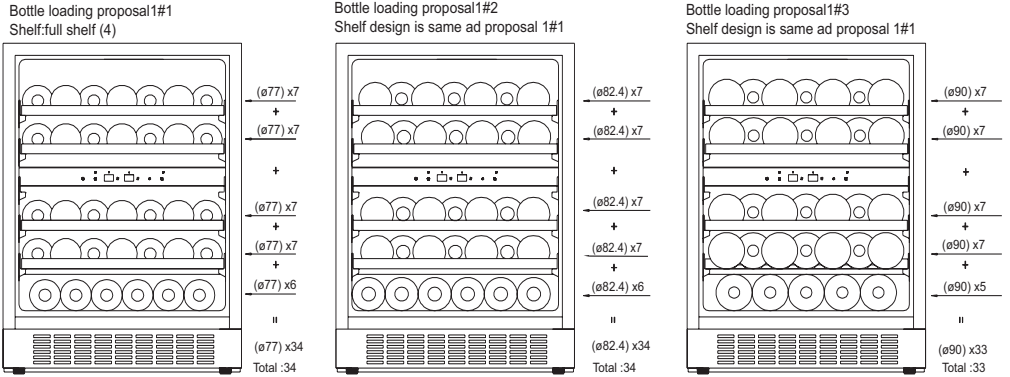
fig. 8

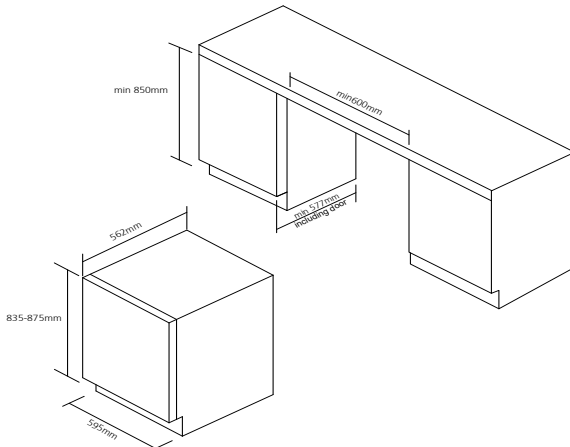
fig. 9

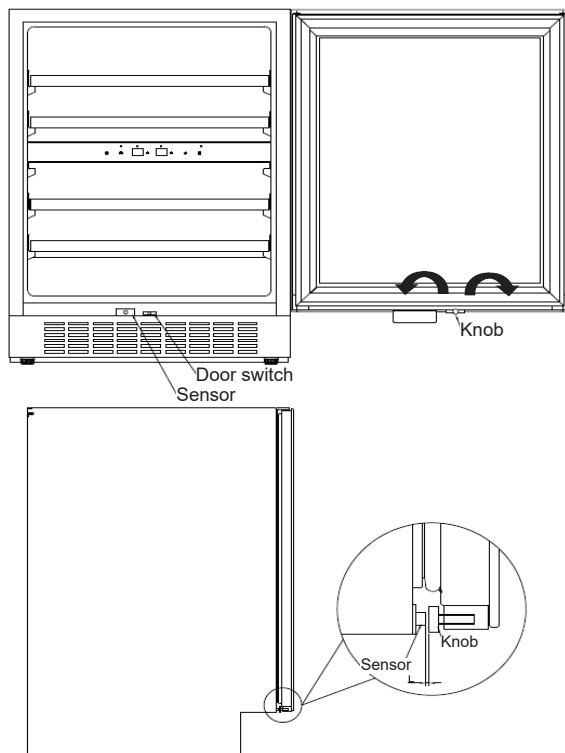
fig. 10

1. The door gasket may be affected by hot expansion which results in less sensitivity. Please follow the instructions to adjust the knob when pushing the door without response
2. When the space between the push-to-open sensor and knob is more than 0.5 mm: Turn the knob to the left, and the knob will extend from door frame. It will extend 0.75 mm when rotated 1 turn.
3. When the space between the push-to-open sensor and knob is less than 0.5 mm: Turn the knob to the right, and the knob will retract to the door frame. It will retract 0.75 mm when rotated 1 turn.

Defrosting, cleaning, and maintenance

Automatic defrosting

The beverage cooler is defrosted automatically. Defrost water runs through a pipe and is collected in a tray above the compressor where the heat generated by the compressor causes it to evaporate.

Cleaning

Before cleaning the appliance, unplug it from the main supply. The cabinet is best cleaned using warm water (max. 65°C) with a little mild detergent. Never use cleaning agents that scour. Use a soft cloth. Rinse with clean water and dry thoroughly. Never use sharp or pointed implements. Cleaning agents may discolour the shelves.

The sealing strip around the door must be cleaned regularly to prevent discolouration and prolong service life. Use clean water. After cleaning the sealing strip, check that it continues to provide a tight seal.

Dust collecting on the condenser underneath the cabinet, on the compressor and in the compressor compartment is best removed using a vacuum cleaner.

Replacing the lighting

Light-emitting diodes designed to last the entire service life of the appliance are used. If it, however, becomes necessary to replace the light source, please call for service.

Fault finding

Fault	Possible cause	Remedy
The appliance is not running	The appliance is switched off. Power failure; the fuse is blown; the appliance is not plugged in correctly.	Press the on/off switch. Check that power is connected. Reset the fuse.
Water collects in the bottom of the cabinet.	The defrost water pipe is blocked.	Clean the defrost water channel and the drain hole on the rear wall of the cabinet.
Vibration or bothersome noise.	The appliance is not level. The appliance is resting against other kitchen elements. Containers or bottles inside the cabinet are rattling against one another.	Level the appliance using a spirit level. Move the appliance away from the kitchen elements or appliances it is in contact with. Move containers and/or bottles apart.
Compressor runs continuously.	High room temperature.	Ensure adequate ventilation.
Beverage cooler is not cold enough	Check the temperature setting. The door is opened too often. The door is not completely closed. The door gasket does not seal properly	Set the desired temperature. Keep the door closed for an extended period of time. Close the door. Align the door.

Warranty, spare parts and service

Warranty disclaimer

Faults and damage caused directly or indirectly by incorrect operation, misuse, insufficient maintenance, incorrect building, installation or mains connection. Fire, accident, lightning, voltage variation or other electrical interference, including defective fuses or faults in mains installations are not covered by the warranty.

Repairs performed by others than approved service centres and any other faults and damage that the manufacturer can substantiate are caused by reasons other than manufacturing or material faults are not covered by the warranty.

Please note that changes to the construction of the appliance or changes to the component equipment of the appliance will invalidate warranty and product liability, and the appliance cannot be used lawfully. The approval stated on rating plate will also be invalidated.

Transport damage discovered by the buyer is primarily a matter to be settled between the buyer and the distributor, i.e. the distributor must ensure that such complaints are resolved to the buyer's satisfaction.

Before calling for technical assistance, please check whether you are able to rectify the fault yourself. (see the fault finding list) If your request for assistance is unwarranted, e.g. if the appliance has failed as a result of a blown fuse or incorrect operation, you will be charged the costs incurred by your call for technical assistance.

Spare parts

When ordering spare parts, please state the model and production numbers of your appliance. This information is given on the rating plate. The rating plate contains various technical information, including type and serial numbers.

VESTFRØST SOLUTIONS Wine cellar WFG48
 MODEL: 0-S710001H8030803

ERIC **CE**

SPECIFICATIONS	
SERIAL NO.:	
CATEGORY:	W
CAPACITY:	8000L
REFRIGERANT:	R135
AMOUNT:	31g
INSULATION:	CFP10000
BLOCK/VALVE:	CFP10000
RATED VOLTAGE:	230/240V
RATED FREQUENCY:	50Hz
RATED CURRENT:	0.6A
NET WEIGHT:	48kg
NET VOLUME:	120L
BOTTLE CAPACITY:	34

FALKEVEJ 12 DK-6705 ESSBJERG, DENMARK
 BUILT-IN OR FREE STANDING INSTALLING

VESTFRØST SOLUTIONS Wine cellar WFG49
 MODEL: 0-S710003H8030803

ERIC **CE**

SPECIFICATIONS	
SERIAL NO.:	
CATEGORY:	W
CAPACITY:	8000L
REFRIGERANT:	R135
AMOUNT:	31g
INSULATION:	CFP10000
BLOCK/VALVE:	CFP10000
RATED VOLTAGE:	230/240V
RATED FREQUENCY:	50Hz
RATED CURRENT:	0.6A
NET WEIGHT:	40kg
NET VOLUME:	120L
BOTTLE CAPACITY:	34

FALKEVEJ 12 DK-6705 ESSBJERG, DENMARK
 BUILT-IN OR FREE STANDING INSTALLING

fig. 11

Note: Always use approved service centres when your appliance is to be repaired or replaced!

Disposal

Information for Users on Collection and Disposal of Old Equipment and used Batteries

These symbols on the products, packaging, and/or accompanying documents mean that used electrical and electronic products and batteries should not be mixed with general household waste. For proper treatment, recovery and recycling of old products and used batteries, please take them to applicable collection points, in accordance with your national legislation and the Directives 2012/19/EU and 2006/66/EC.

By disposing of these products and batteries correctly, you will help to save valuable resources and prevent any potential negative effects on human health and the environment which could otherwise arise from inappropriate waste handling.

For more information about collection and recycling of old products and batteries, please contact your local municipality, your waste disposal service or the point of sale where you purchased the items.

Penalties may be applicable for incorrect disposal of this waste, in accordance with national legislation.

For business users in the European Union.

If you wish to discard electrical and electronic equipment, please contact your dealer or supplier for further information.

Information on Disposal in other Countries outside the European Union

These symbols are only valid in the European Union. If you wish to discard this product, please contact your local authorities or dealer and ask for the correct method of disposal.

Note for the battery symbol:

This symbol might be used in combination with a chemical symbol. In this case it complies with the requirement set by the Directive for the chemical involved.

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A/S VESTFROST
A/S Vestfrost